

Zeta Plus™ VR Series

Depth Filter Cartridges & Capsules

Application

The removal and/or inactivation of contaminating viruses from biotherapeutics is a requisite for ensuring final product safety. Zeta Plus VR Series cartridge depth filters remove significant levels of viruses from biological fluids. They provide validatable viral titer reduction, high flow rates, scalability, economy, disposability and ease-of-use in the biological manufacturing environment. The Zeta Plus VR Series includes specific filter media recommendations for virus removal from blood plasma proteins and bioprocess-derived cell culture fluids.

Virus Removal Applications for Zeta Plus VR Media:

- Processing of human plasma-derived proteins
- Recombinant proteins from cell culture and fermentation
- Monoclonal antibodies
- Serum and other raw materials derived from animal sources

Zeta Plus depth filtration offers an attractive and economical alternative to other available technologies for reduction of viral burden. Zeta Plus VR Series depth filters can be used as a bolt-on processing step to existing in-process viral clearance methodologies, providing an added measure of end-product safety.



Zeta Plus VR Series For
Enhanced Log Reduction
of Viruses From Biological
Fluids

Features & Benefits

Combined ion exchange & physical entrapment virus removal mechanisms.

- Validatable viral log reduction.
- Provides added level of viral safety.

Economical “bolt-on” separation technology as adjunct or pre-filter to other viral clearance steps.

- Provides added level of validatable viral clearance.
- Can replace or augment chromatography columns or function as prefilter to final virus removal membrane filter.

Quality Control tested for media ion exchange capacity.

- Assurance of filter media consistency & performance.

Self-contained, single-use disposable capsules & cartridge modules

- Single-use, disposable design eliminates validation concerns associated with reusable clearance options (chromatography column sanitization & regeneration).

Range of cartridge sizes, sanitary design cartridge housings & disposable capsules.

- Facilitates scaleup from small volume feasibility trials & validation studies to production scale operations.



Zeta Plus VR Series filter media is available in a wide range of cartridge sizes including 8-, 12- and 16-inch diameter filter cartridges for varying flow rates and batch volumes. 3M Purification Inc.'s sanitary design ZPB Zeta Plus filter housings are designed for pharmaceutical and biological manufacturing environments. For more information on Zeta Plus sanitary filter housings, request 3M Purification Inc. literature numbers 70-0201-8857-2 and 70-0201-8762-4.

Zeta Plus VR Series filter media is a family of cellulosic depth filtration media designed to retain contaminants by both physical entrapment and ion exchange adsorption. It is composed of high-area process filter aids embedded in a cellulose fiber depth matrix. During the manufacturing process, a cationic charge modifier is chemically bound to the matrix components, forming a permanent, interconnected, rigid depth filter with positively-charged electrokinetic capture sites. The resulting porous depth filter structure is a tortuous network of adsorptive flow channels capable of retaining contaminating viruses through a combination of anion exchange adsorption and mechanical entrapment. Cumulative viral titer reduction may also be attainable through a staged two-step Zeta Plus system.

Zeta Plus depth filtration is a key component in biopharmaceutical production schemes for clarifying cell culture fluids (cell separation) and is utilized world wide in the fractionation of human and animal plasma proteins.

Zeta Plus VR Series Filter Selection

A range of cartridge sizes is available for varying batch size, so scale-up is easy and predictable from lab-scale, to pilot plant, to full production scale. Three different cartridge sizes (8, 12 and 16-inch diameter) make scale-up straight forward and predictable. Zeta Plus cartridges are used in a flow-through mode, enabling the process stream to be passed through the filter cartridge while viruses are retained within the depth filter matrix. Small, low area disposable BC capsules are available for laboratory scale and process development work and are ideal for scaled down viral validation studies.

Zeta Plus VR Series Disposable Capsules, Filter Cartridges, and Filter Sheets—Configured to Suit Your Every Need

Zeta Plus VR Series filter products are highly scaleable and can be used for applications ranging from milliliter lab volumes to production scale cGMP operations. Filter media is available in the following configurations:

Zeta Plus VR Series Filter Sheets: VR filter media is available in a range of filter sheet sizes to fit plate-and-frame filter presses (contact 3M Purification Inc. for specific sheet sizes and part numbers).

Zeta Plus VR Series BC Disposable Filter Capsules: VR filter media is available in three different sizes of BC disposable capsules for bench scale and pilot scale filtration.

Zeta Plus VR Series Filter Cartridges and Sanitary Housings: VR filter media is available in 8-inch, 12-inch and 16-inch diameter, easy-to-use, disposable cartridges. A range of sanitary design Zeta Plus filter housings is available for VR Series cartridges.


Table 1: Selection of Zeta Plus VR Series Filter Media

| Filter Media Choice | Recommended For: | | | | | |
|---------------------|------------------|-----------------------|-------------------------------|------------------------------------|--|-------------------------------------|
| | Plasma Proteins | Cell Culture Proteins | Nominal Retention Rating (µm) | 20°C Water Flow @ 5 psid (gpm/ft²) | Autoclavable or In-situ Steam Sterilizable | [AI3+] in 20°C water @ 5L/ft² (ppb) |
| VR 01 | • | | 0.8 | 1.2 | No | 40 |
| VR 02 | • | • | 0.8 | 1.9 | Yes | 20 |
| VR 03 | | • | 0.8 | 1.2 | Yes | 40 |
| VR 04 | • | | 0.45 | 0.9 | No | 20 |
| VR 05 | | • | 0.2 | 0.4 | Yes | 40 |
| VR 06 | • | • | 0.2 | 0.6 | Yes | 20 |
| VR 07 | • | • | 0.2 | 0.4 | Yes | 60 |

Recommended Operating Parameters

| | |
|-------------------------------|---|
| Nominal Retention Ratings | 0.2 to 0.8 micron |
| Maximum Operating Pressure | VR Cartridges: 35 psi maximum cartridge pressure drop. BC25 Capsules: 40 psig max. inlet pressure, 35 psi max. media pressure drop. BC1000/2000 Series Capsules: 80 psig at 25° C, 40psig at 60° C, max. inlet pressure; 35 psi max. cartridge pressure drop. |
| Maximum Operating Temperature | Zeta Plus VR 02, 03, 05, 06, & 07 Cartridges: 82° C. Zeta Plus VR 01 & 04 Cartridges: 60° C. BC25 Capsules: 40° C. BC1000 / 2000 Capsules: 60° C. |
| Recommended Pre-use Rinse | 1.3 gal/ft² at 0.5 gpm/ ft², 50 L/m² at 20 L/m²/min. |
| Sterilization Parameters | All media grades can be autoclaved or <i>in-situ</i> steam sterilized, except VR 01 & VR 04 (refer to Table 1). Zeta Plus VR Series Cartridges: In-situ steam sterilized or autoclaved, for 30 minutes at 121°C (1 cycle). BC25: Autoclave for 30 minutes at 121°C (1 cycle). BC1000/2000: Autoclave for 30 minutes at 121°C (up to 3 cycles). |
| Recommended Flow Rates | BC25: 3 - 30 ml/min. BC1000/2000 & VR Cartridges: 0.065 gpm/ft², 2.5 L/m²/min. |

Cartridge Construction

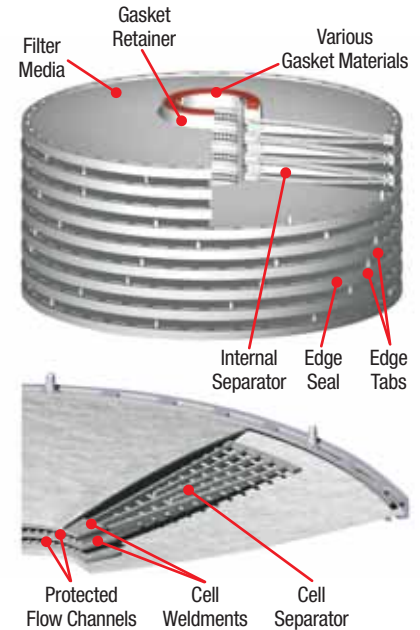

Viral Reduction Validation

Many factors influence the choice of process steps to be studied when performing viral clearance evaluation and validation. Viral clearance validation studies are performed with a select panel of model viruses. Different model viruses are chosen to validate blood and plasma products, and separately, cell-culture derived products.

Virus removal efficiency may be dependent on fluid and processing conditions and is best determined during qualification and validation studies using suitable virus model systems in a controlled laboratory study. Contact 3M Purification Inc. Scientific Applications Support Services (SASS) for further technical advice concerning validation of viral reduction with Zeta Plus VR Series depth filters.

Data is available which confirms the ability of Zeta Plus VR Series depth filter media to function as an effective and validatable viral log reduction filter. Table 2 illustrates results for a reported viral clearance study utilizing Zeta Plus VR depth filters.

Quality Control of Zeta Plus VR Series Filters

3M Purification Inc. applies rigorous Quality Control testing and standards both during in-process manufacturing and during final lot release. This testing ensures consistent filter media performance in critical virus removal applications. A Certificate of Quality is provided with each Zeta Plus VR Series filter cartridge and disposable BC capsule.

Each grade of Zeta Plus VR Series filter media is Quality Control tested on a lot release basis for the presence and magnitude of positive charge.

Table 2: Viral Clearance In An Immunoglobulin Production Process
Cumulative Virus Titer Reduction (Log₁₀)

| Process Step | BVD | EMC | HIV | PPV | PRV |
|-----------------------------------|-----------------|------------|-----------------|------------|-----------------|
| Solvent detergent | >4.3 | — | >5.3 | — | >7.3 |
| Supernatant III | 1.4 | 4.3 | 6.1 | 4.7 | 3.8 |
| Zeta Plus VR 03 Depth Filtration | 4.8 | 4.5 | 4.7 | 3.7 | 5.4 |
| Total Cumulative Reduction | >10.5 | 8.8 | >16.1 | 8.4 | >16.6 |

Source: D. Revie, *Novel Validation Approaches to Obtain Maximum Viral Clearance from an Immunoglobulin Production Process*, IBC 2nd International Symposium on Viral Clearance, Philadelphia, PA, June, 1998.

| Effective Filtration Area | |
|---|--------------------|
| BC25 Capsule | 3.9 in² (25 cm²) |
| BC1000 Capsule | 0.7 ft² (650 cm²) |
| BC2000 Capsule | 1.4 ft² (1300 cm²) |
| 45109 (8-inch diameter cartridge, 8-cell) | 2.8 ft² (0.26 m²) |
| 45167 (8-inch diameter cartridge, 7-cell, o-ring plug-in) | 2.5 ft² (0.23 m²) |
| Z8FA2NPX2 (8-inch diameter cartridge, 2-cell plug-in) | 0.7 ft² (650 cm²) |
| Z8FA4NPX2 (8-inch diameter cartridge, 4-cell plug-in) | 1.4 ft² (1300 cm²) |
| 45264 (12-inch diameter cartridge, 7-cell bodyfeed) | 7.2 ft² (0.67 m²) |
| 45245 (12-inch diameter cartridge, 16-cell) | 16.4 ft² (1.5 m²) |
| Z16P (16-inch, 14-cell, netted) | 34.7 ft² (3.2 m²) |

Zeta Plus™ VR Series Ordering Guide

8" Diameter Cartridges

| Catalog Number | Configuration | VR Series |
|--|--|---|
| 45109 (8" 8-cell) | 11 - Nitrile 13 - Fluorocarbon 14 - EPR 22 - Silicone 23 - Expanded PTFE | VR01 VR05 VR02 VR06 VR03 VR07 VR04 |
| 45167 (8" 7-cell, O-ring plug-in) | 01 - Nitrile 02 - EPR 03 - Fluorocarbon 04 - Silicone | VR01 VR05 VR02 VR06 VR03 VR07 VR04 |
| Z8FA2NP (8" 2-cell, plug-in) Z8FA4NP (8" 4-cell, plug-in) | 01 - Nitrile 02 - EPR 03 - Fluorocarbon 04 - Silicone | VR01 VR05 VR02 VR06 VR03 VR07 VR04 |
| | Pkg Code 2 | |

16" Diameter Cartridges

| Catalog Number | Gasket Material | VR Series |
|--------------------|--|---|
| Z16P (16" 14-cell) | A - Silicone B - Fluorocarbon C - EPR D - Nitrile | VR01 VR05 VR02 VR06 VR03 VR07 VR04 |

BC Capsules

| Catalog Number | VR Series |
|--|---|
| BC0025L (Luer) BC0025S (Sanitary) BC1000A (Single pack) BC1000B (3 pack) BC2000A (Single pack) BC2000B (3 pack) | VR01 VR05 VR02 VR06 VR03 VR07 VR04 |

12" Diameter Cartridges

| Catalog Number | Geometric Variation | Gasket Material | VR Series |
|---|--------------------------------|--|---|
| 45264 (12" 7-cell, Bodyfeed) 45245 (12" 16-cell) | 01 - Standard Polypropylene | A - Silicone B - Fluorocarbon C - EPR D - Nitrile | VR01 VR05 VR02 VR06 VR03 VR07 VR04 |

Important Notice: The information described in this literature is accurate to the best of our knowledge. A variety of factors, however, can affect the performance of the Product(s) in a particular application, some of which are uniquely within your knowledge and control. **INFORMATION IS SUPPLIED UPON THE CONDITION THAT THE PERSONS RECEIVING THE SAME WILL MAKE THEIR OWN DETERMINATION AS TO ITS SUITABILITY FOR THEIR USE. IN NO EVENT WILL 3M PURIFICATION INC. BE RESPONSIBLE FOR DAMAGES OF ANY NATURE WHATSOEVER RESULTING FROM THE USE OF OR RELIANCE UPON INFORMATION.**

It is your responsibility to determine if additional testing or information is required and if this product is fit for a particular purpose and suitable in your specific application.

3M PURIFICATION INC. MAKES NO REPRESENTATIONS OR WARRANTIES, EITHER EXPRESS OR IMPLIED INCLUDING WITHOUT LIMITATION ANY WARRANTIES OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE OR OF ANY OTHER NATURE HEREUNDER WITH RESPECT TO INFORMATION OR THE PRODUCT TO WHICH INFORMATION REFERS

Limitation of Liability: 3M Purification Inc. will not be liable for any loss or damage arising from the use of the Product(s), whether direct, indirect, special, incidental, or consequential, regardless of the legal theory asserted, including warranty, contract, negligence or strict liability. Some states do not allow the exclusion or limitation of incidental or consequential damages, so the above limitation may not apply to you.

Your Local 3M Purification Inc. Distributor:



3M Purification Inc.

400 Research Parkway
Meriden, CT 06450
U.S.A.
(800) 243-6894
(203) 237-5541
Fax (203) 630-4530
www.3mpurification.com

3M is a trademark of 3M Company.
Zeta Plus is a trademark of 3M Company used under license.
© 2011 3M Company. All rights reserved.
70-0201-8875-4
REV 0411pb